



DC1500 CHLORINE



Chlorine Colorimeter Kit • DPD Liquid

Code 3240-LI

Firmware Version 2.10 or greater

CONTENTS

Chlorine & Bromine DPD Liquid Kit	
■ Kit Contents.....	5
■ Accessories	5
Test Method Specifications	
■ Application.....	6
■ Range.....	6
■ Method	6
■ Handling & Preservation	6
■ Interferences.....	6
■ Calibration	6
Analysis	
■ Chlorine	7
■ Bromine.....	10
Calibration	
■ Chlorine Standards	12
■ Chlorine Calibration Procedure	12
Set Up	
■ Setting the Clock	15
■ Logging Data	16
■ Factory Setup.....	17
■ Setting Power Save	17
■ Setting Backlight Time	18
■ Selecting A Language	19
■ Looping Menus.....	20
Computer Connection	
■ Output.....	21
■ Computer Connection.....	21
Battery	
■ Battery/AC Operation.....	21
■ Battery Replacement.....	22
Maintenance	
■ Cleaning	22
■ Repairs	22
■ Meter Disposal	23
General Operating Information	
■ Overview.....	24
■ General Operating Information	24
■ The Keypad	24
■ The Display and Menus	25

■ Tubes and Chambers.....	27
■ Sample Dilution Techniques	27
General information	
■ Packaging and Delivery	28
■ General Precautions.....	28
■ Safety Precautions.....	28
■ Limits of Liability	29
■ Specifications	29
■ CE	30
■ IP67.....	30
■ Warranty	30
■ Register Your Meter.....	30
Trouble Shooting Guide	
■ Error Messages	31
■ Calibration	31
■ Stray Light	31
■ Troubleshooting Guide.....	32

DC1500 Chlorine & Bromine DPD Liquid Kit

■ Kit Contents

QUANTITY	CONTENTS	CODE
30 mL	DPD 1A Free Chlorine Reagent	P-6740-G
30 mL	*DPD 1B Free Chlorine Reagent	*P-6741-G
30 mL	*DPD 3 Total Chlorine Reagent	*P-6743-G
1	Colorimeter Tubes, w/caps, set of 6	0290-6
1	Water Sample Collecting Bottle	0688
1	1500 Colorimeter for Chlorine DPD	27926-CL
1	USB Wall Adapter	1721
1	USB Cable	1720
1	DC1500-CL Chlorine Colorimeter, DPD Liquid Manual	3240-LI-MN-CL
1	DC1500-CL Chlorine Colorimeter, DPD Liquid Quick Start Guide	3240-LI-QG-CL

***WARNING:** Reagents marked with an * are considered to be potential health hazards. To view or print a Safety Data Sheet [SDS] for these reagents go to [www.itm.com](#). Search the four digit reagent code number listed on the reagent label, in the contents list or in the test procedures. Omit any letter that follows or precedes the four digit code number. For example if code is 4450WT-H, search 4450. To obtain a printed copy, contact LaMotte by e-mail, phone or fax.

To order individual reagents or test kit components, use the specified code number.

■ Accessories

DESCRIPTION	CODE
Test Tubes, with Caps	0290-6
Replacement Chamber	3-0038
USB Cable	1720
USB Wall Adapter	1721
Car Charger	5-0132

TEST METHODS SPECIFICATIONS

■ APPLICATION

Drinking water supplies and distribution systems, swimming pool and spas, sewage and chlorinated waste waters, process waters and sanitizing solutions.

■ RANGE

0 to 4.0 ppm Chlorine [may be extended by dilution]

0 to 7.0 ppm Bromine [may be extended by dilution]

■ METHOD

In the absence of Iodide, Free Available Chlorine reacts instantly with the buffered diethyl-p-phenylenediamine indicator (DPD) to produce a pink-red color in proportion to the amount of chlorine present. Subsequent addition of potassium iodide produces a rapid color response from the combined forms of chlorine [chloramines]. In buffered samples, bromine reacts with diethyl-p-phenylene diamine (DPD) to produce a pink-red color in proportion to the concentration of bromine present.

■ HANDLING & PRESERVATION

Chlorine and bromine in aqueous solutions, particularly weak solutions, are not stable. Exposure to sunlight or agitation will accelerate the reduction of chlorine and bromine. For chlorine samples, fill sample containers to the top and cap tightly. Analyze samples as soon as possible after collection. Samples to be analyzed for bromine cannot be preserved or stored.

■ INTERFERENCES

The only interfering substance likely to be encountered is oxidized manganese. The extent of this interference can be determined by treating a sample with sodium arsenite to destroy the chlorine or bromine present, so that the amount of interference can be estimated. Chlorine, bromine, and iodine will give a positive result and should be considered an interference unless the test is being performed specifically for that parameter.













■ CALIBRATION

The colorimeter is precalibrated. In order to comply with NPDWR or NPDES reporting regulations, the calibration should be checked periodically by using a set of reference standards including a 0 mg/L blank and 0.3, 1.0, and 3.5 mg/L chlorine. To prepare these standards, a LaMotte 1000 mg/L standard chlorine equivalent solution [Code 3858] is available. Consult with your local regulatory agency to determine standardization frequency.


ANALYSIS - CHLORINE

PROCEDURE – FREE CHLORINE


Chlorine

<p>1. Fill the Water Sample Collecting Bottle (0688) with sample water.</p>																			
<p>2. Press and hold  until colorimeter turns on.</p>	<table border="1"> <tr><th colspan="3">Main Menu</th></tr> <tr><td colspan="3">Testing Menu</td></tr> <tr><td colspan="3">Editing Menu</td></tr> <tr><td>12:00:00</td><td>001/500</td><td></td></tr> </table>	Main Menu			Testing Menu			Editing Menu			12:00:00	001/500							
Main Menu																			
Testing Menu																			
Editing Menu																			
12:00:00	001/500																		
<p>3. Press ENTER to select Testing Menu.</p>	<table border="1"> <tr><th colspan="3">All Tests</th></tr> <tr><td colspan="3">015 Bromine L DPD</td></tr> <tr><td colspan="3">016 Bromine T DPD</td></tr> <tr><td colspan="3">024 Chlorine L DPD</td></tr> <tr><td colspan="3">025 Chlorine T DPD</td></tr> <tr><td>12:00:00</td><td>001/500</td><td></td></tr> </table>	All Tests			015 Bromine L DPD			016 Bromine T DPD			024 Chlorine L DPD			025 Chlorine T DPD			12:00:00	001/500	
All Tests																			
015 Bromine L DPD																			
016 Bromine T DPD																			
024 Chlorine L DPD																			
025 Chlorine T DPD																			
12:00:00	001/500																		
<p>4. Scroll to and select 024 Chlorine L DPD from menu.</p>	<table border="1"> <tr><th colspan="3">024 Chlorine L DPD</th></tr> <tr><td colspan="3">[blacked out]</td></tr> <tr><td colspan="3">Scan Blank</td></tr> <tr><td colspan="3">Scan Sample</td></tr> <tr><td>12:00:00</td><td>001/500</td><td></td></tr> </table>	024 Chlorine L DPD			[blacked out]			Scan Blank			Scan Sample			12:00:00	001/500				
024 Chlorine L DPD																			
[blacked out]																			
Scan Blank																			
Scan Sample																			
12:00:00	001/500																		
<p>5. Rinse a clean tube (0290) with sample water. Fill to the 10 mL line with sample. Cap and wipe dry.</p>																			

6. Insert tube into chamber, close lid and select **Scan Blank**.

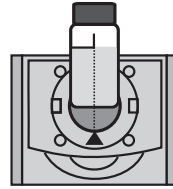

7. Remove tube from colorimeter. Add 5 drops of DPD 1A Free Chlorine Reagent (P-6740) and add 5 drops of *DPD 1B Free Chlorine Reagent (P-6741), Cap and invert to mix. The solution will be pink if chlorine is present. Read within 30 seconds.

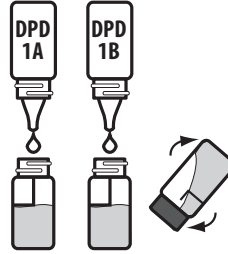

8. Immediately insert tube into chamber. Close lid. Select **Scan Sample**. Record result as Free Chlorine.

024 Chlorine L DPD

0.99 ppm

Scan Blank

Scan Sample


12:00:00

001/500




PROCEDURE – COMBINED CHLORINE

9. Remove tube from colorimeter. Add 5 drops of *DPD 3 Total Chlorine Reagent (P-6743). Cap and invert to mix. An increase in color represents Combined Chlorine.

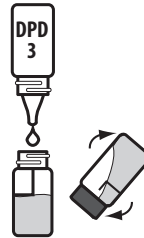



NOTE: For wastewater samples, *Standard Methods for the Examination of Water and Wastewater* recommends waiting 2 minutes for full color development when testing total chlorine.

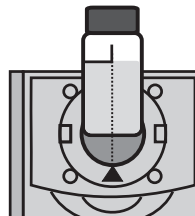10. Insert tube into chamber, close lid and select **Scan Sample**. Record result as Total Chlorine (Monochloramine, Dichloramine, and Nitrogen Trichloride) present in the water sample.



11. Subtract free chlorine reading from total chlorine reading to obtain concentration of combined chlorine.

12. Press  to turn off the colorimeter or press  to exit to a previous menu or make another menu selection.



















Levels of chlorine about 4 mg/L will cause a bleaching effect on the DPD indicator, and may give a false indication that no chlorine is present. If it is possible that the chlorine concentration is greater than 4 mg/L, perform test on a diluted sample and multiply the result by appropriate dilution factor.


CAUTION: DO NOT leave reacted samples in test tubes. Discard sample and thoroughly rinse tubes. If allowed to remain, DPD will stain tubes, significantly impairing the operation of the colorimeter. If necessary, acid wash, and vigorously clean glassware with test tube brush and detergent.


NOTE: For the most accurate results, samples over 4 ppm chlorine should be diluted with chlorine demand free water and re-tested.




NOTE: The meter will remember the last scanned blank reading. It is not necessary to scan a blank each time the test is performed. To use the previous blank reading, instead of scanning a new one, scroll to Scan Sample and proceed. For the most accurate results, the meter should be blanked before each test and the same tube should be used for the blank and the reacted sample.

ANALYSIS - BROMINE PROCEDURE
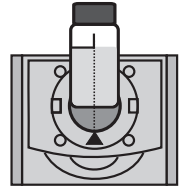
<p>1. Fill the Water Sample Collecting Bottle (0688) with sample water.</p>																			
<p>2. Press and hold  until colorimeter turns on.</p>	<table border="1"> <tr> <td colspan="3">Main Menu</td> </tr> <tr> <td colspan="3">Testing Menu</td> </tr> <tr> <td colspan="3">Editing Menu</td> </tr> <tr> <td>12:00:00</td> <td>001/500</td> <td></td> </tr> </table>	Main Menu			Testing Menu			Editing Menu			12:00:00	001/500							
Main Menu																			
Testing Menu																			
Editing Menu																			
12:00:00	001/500																		
<p>3. Press ENTER to select Testing Menu.</p>	<table border="1"> <tr> <td colspan="3">All Tests</td> </tr> <tr> <td colspan="3">015 Bromine L DPD</td> </tr> <tr> <td colspan="3">016 Bromine T DPD</td> </tr> <tr> <td colspan="3">024 Chlorine L DPD</td> </tr> <tr> <td colspan="3">025 Chlorine T DPD</td> </tr> <tr> <td>12:00:00</td> <td>001/500</td> <td></td> </tr> </table>	All Tests			015 Bromine L DPD			016 Bromine T DPD			024 Chlorine L DPD			025 Chlorine T DPD			12:00:00	001/500	
All Tests																			
015 Bromine L DPD																			
016 Bromine T DPD																			
024 Chlorine L DPD																			
025 Chlorine T DPD																			
12:00:00	001/500																		
<p>4. Scroll to and select 015 Bromine L DPD from menu.</p>	<table border="1"> <tr> <td colspan="3">015 Bromine L DPD</td> </tr> <tr> <td colspan="3"></td> </tr> <tr> <td colspan="3">Scan Blank</td> </tr> <tr> <td colspan="3">Scan Sample </td> </tr> <tr> <td>12:00:00</td> <td>001/500</td> <td></td> </tr> </table>	015 Bromine L DPD						Scan Blank			Scan Sample 			12:00:00	001/500				
015 Bromine L DPD																			
																			
Scan Blank																			
Scan Sample 																			
12:00:00	001/500																		
<p>5. Rinse a clean tube (0290) with sample water. Fill to the 10 mL line with sample. Cap and wipe dry.</p>																			
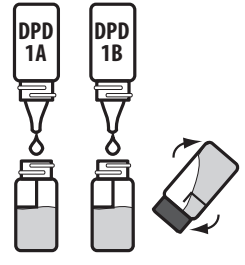
<p>6. Insert tube into chamber, close lid and select Scan Blank.</p>	
---	--

<p>7. Remove tube from colorimeter. Add 5 drops of DPD 1A Free Chlorine Reagent (P-6740) and add 5 drops of *DPD 1B Free Chlorine Reagent (P-6741), Cap and invert to mix. The solution will be pink if bromine is present. Read within 30 seconds.</p>	
---	--

<p>8. Immediately insert tube into chamber. Close lid. Select Scan Sample. Record result as Bromine.</p>	<table border="1"> <tr> <td colspan="2" style="text-align: center;">015 Bromine L DPD</td> </tr> <tr> <td colspan="2" style="text-align: center; font-size: 2em;">0.99 ppm</td> </tr> <tr> <td colspan="2" style="text-align: center;">Scan BLank</td> </tr> <tr> <td colspan="2" style="text-align: center;">Scan Sample</td> </tr> <tr> <td style="text-align: center;">12:00:00</td> <td style="text-align: center;">001/500 </td> </tr> </table>	015 Bromine L DPD		0.99 ppm		Scan BLank		Scan Sample		12:00:00	001/500 
015 Bromine L DPD											
0.99 ppm											
Scan BLank											
Scan Sample											
12:00:00	001/500 										

NOTE: For the most accurate results, samples over 7 ppm bromine should be diluted with chlorine demand free water and re-tested.

NOTE: The meter will remember the last scanned blank reading. It is not necessary to scan a blank each time the test is performed. To use the previous blank reading, instead of scanning a new one, scroll to Scan Sample and proceed. For the most accurate results, the meter should be blanked before each test and the same tube should be used for the blank and the reacted sample.

CALIBRATION


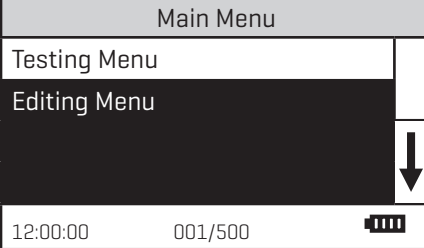
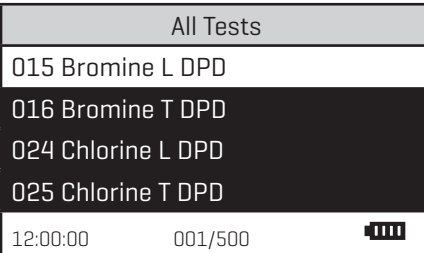
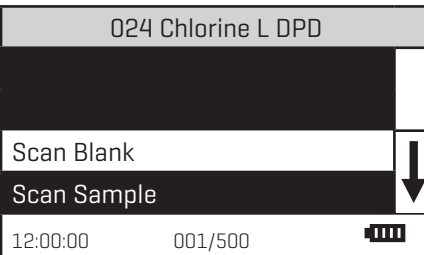
■ CHLORINE STANDARDS


The meter should be calibrated with free chlorine standards. The calibration should be done with a distilled or deionized water blank and one chlorine standard of known concentration. The concentration of the calibration standard should be similar to the expected concentration of the sample that will be tested.


To perform a bromine calibration, prepare a bromine with a concentration that is equivalent to a chlorine dioxide standard that is within the range of the chlorine dioxide test. (1 ppm chlorine = 1.9 ppm bromine). Follow the procedure below but select the bromine test.


Chlorine Secondary Standards [Code 4140-03] are available to verify the performance of the meter.




CHLORINE CALIBRATION PROCEDURE



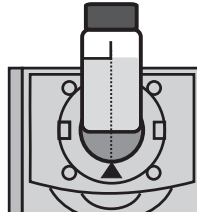
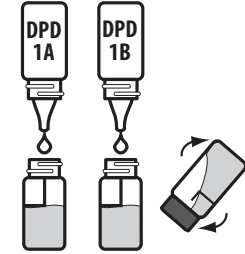
<p>1. Press and briefly hold  to turn the meter on. The LaMotte logo screen will appear for about 3 seconds and the Main Menu will appear.</p>	 <p>The screenshot shows the 'Main Menu' with options 'Testing Menu' and 'Editing Menu'. A downward arrow is on the right side. At the bottom, it displays '12:00:00', '001/500', and a battery level icon.</p>
<p>2. Press ENTER to select Testing Menu.</p>	 <p>The screenshot shows the 'All Tests' menu with options: '015 Bromine L DPD', '016 Bromine T DPD', '024 Chlorine L DPD', and '025 Chlorine T DPD'. A downward arrow is on the right side. At the bottom, it displays '12:00:00', '001/500', and a battery level icon.</p>
<p>3. Scroll to and select 024 Chlorine L DPD from menu.</p>	 <p>The screenshot shows the '024 Chlorine L DPD' menu with options 'Scan Blank' and 'Scan Sample'. A downward arrow is on the right side. At the bottom, it displays '12:00:00', '001/500', and a battery level icon.</p>





<p>4. Rinse a clean tube (0290) three times with the chlorine standard. Fill to the 10 mL line with the chlorine standard. Cap the tube. Dry the tube with a lint-free cloth.</p>	
---	--

<p>5. Insert tube into chamber, close lid and select Scan Blank.</p>	
---	--

<p>6. Remove tube from colorimeter. Add 5 drops of DPD 1A Free Chlorine Reagent (P-6740) and add 5 drops of *DPD 1B Free Chlorine Reagent (P-6741), Cap and invert to mix. The solution will be pink if chlorine is present. Read within 30 seconds.</p>	
--	---

<p>7. Immediately insert tube into chamber. Close lid. Select Scan Sample.</p>	<table border="1"> <tr> <td colspan="2">024 Chlorine L DPD</td> </tr> <tr> <td colspan="2">0.99 ppm</td> </tr> <tr> <td>Scan BLank</td> <td>↑</td> </tr> <tr> <td>Scan Sample</td> <td>↓</td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </table>	024 Chlorine L DPD		0.99 ppm		Scan BLank	↑	Scan Sample	↓	12:00:00	001/500 
024 Chlorine L DPD											
0.99 ppm											
Scan BLank	↑										
Scan Sample	↓										
12:00:00	001/500 										























<p>8. Press  to scroll to Calibrate.</p>	<table border="1"> <tr> <td colspan="2">024 Chlorine L DPD</td> </tr> <tr> <td colspan="2">0.99 ppm</td> </tr> <tr> <td>%T/Abs</td> <td>↑</td> </tr> <tr> <td>Calibrate</td> <td>↓</td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </table>	024 Chlorine L DPD		0.99 ppm		%T/Abs	↑	Calibrate	↓	12:00:00	001/500 
024 Chlorine L DPD											
0.99 ppm											
%T/Abs	↑										
Calibrate	↓										
12:00:00	001/500 										

<p>9. Press ENTER to select Calibrate. A reverse font (light background with dark characters) will appear to indicate that the reading can be adjusted.</p>	<p>024 Chlorine L DPD</p> <p>0.99 ppm</p> <p>^, v=Edit, ENTER=Save ^ +ENTER=Default</p> <p>12:00:00 001/500 </p>	
	<p>10. Press ^ or v to adjust the value shown to the concentration of the prepared standard, 1.00 in this example. NOTE: A maximum adjustment of 25% is possible.</p>	<p>024 Chlorine L DPD</p> <p>1.00 ppm</p> <p>^, v=Edit, ENTER=Save ^ +ENTER=Default</p> <p>12:00:00 001/500 </p>
	<p>11. Press ENTER to save the value.</p> <p>To leave the Calibration procedure without saving the adjustment, press EXIT.</p> <p>Press ^ and ENTER at any time to return to the default value.</p> <p>The calibration has now been standardized and can be used for testing. Scroll to Scan Blank and begin testing.</p>	<p>024 Chlorine L DPD</p> <p>1.00 ppm </p> <p>%T/Abs</p> <p>Calibrate</p> <p>12:00:00 001/500 </p>

SET UP



































































■ SETTING THE CLOCK

Setting the clock allows the correct time and date stamp to be stored with each reading in the data logger.

<p>1. From the Editing Menu, press  or  to scroll to Set Clock.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Set Clock</td> <td rowspan="4" style="text-align: center; vertical-align: middle;">   </td> </tr> <tr> <td>Logging</td> </tr> <tr> <td>Factory Setup</td> </tr> <tr> <td>Set Power Save</td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Editing Menu		Set Clock	 	Logging	Factory Setup	Set Power Save	12:00:00 001/500 				
Editing Menu													
Set Clock	 												
Logging													
Factory Setup													
Set Power Save													
12:00:00 001/500 													
<p>2. Press  to select Set Clock. The year is displayed. Press  or  to scroll to the appropriate character. Press  to select the character. The month, day, hour, format hour, minute, second, AM/PM will be displayed. Repeat for each.</p>	<table border="1"> <thead> <tr> <th colspan="2">Set Time</th> </tr> </thead> <tbody> <tr> <td>Year: 20<u>00</u></td> <td></td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Set Time		Year: 20 <u>00</u>		12:00:00 001/500 							
Set Time													
Year: 20 <u>00</u>													
12:00:00 001/500 													
<p>3. Press  to select the final character. The time and date will be saved and the meter will return to the Editing Menu.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Set Clock</td> <td></td> </tr> <tr> <td>Logging</td> <td></td> </tr> <tr> <td>Factory Setup</td> <td></td> </tr> <tr> <td>Set Power Save</td> <td></td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Editing Menu		Set Clock		Logging		Factory Setup		Set Power Save		12:00:00 001/500 	
Editing Menu													
Set Clock													
Logging													
Factory Setup													
Set Power Save													
12:00:00 001/500 													

■ LOGGING DATA

The default setting for the data logger is enabled. The meter will log the last 500 data points. The counter in the center bottom of the display will show how many data points have been logged. The display will show 500+ when the data logger has exceeded 500 points and the data points are being overwritten.































<p>1. From the Editing Menu, press  or  to scroll to Logging.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Set Clock</td> <td></td> </tr> <tr> <td>Logging</td> <td></td> </tr> <tr> <td>Factory Setup</td> <td></td> </tr> <tr> <td>Set Power Save</td> <td></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Editing Menu		Set Clock		Logging		Factory Setup		Set Power Save		12:00:00	001/500 
Editing Menu													
Set Clock													
Logging													
Factory Setup													
Set Power Save													
12:00:00	001/500 												
<p>2. Press  to select Logging.</p>	<table border="1"> <thead> <tr> <th colspan="2">Logging</th> </tr> </thead> <tbody> <tr> <td>Display Test Log</td> <td></td> </tr> <tr> <td>Logging Enabled</td> <td></td> </tr> <tr> <td>Logging Disabled</td> <td></td> </tr> <tr> <td>Erase Log</td> <td></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Logging		Display Test Log		Logging Enabled		Logging Disabled		Erase Log		12:00:00	001/500 
Logging													
Display Test Log													
Logging Enabled													
Logging Disabled													
Erase Log													
12:00:00	001/500 												
<p>3. Press  or  to scroll to desired function.</p>	<table border="1"> <thead> <tr> <th colspan="2">Logging</th> </tr> </thead> <tbody> <tr> <td>Display Test Log</td> <td></td> </tr> <tr> <td>Logging Enabled</td> <td></td> </tr> <tr> <td>Logging Disabled</td> <td></td> </tr> <tr> <td>Erase Log</td> <td></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Logging		Display Test Log		Logging Enabled		Logging Disabled		Erase Log		12:00:00	001/500 
Logging													
Display Test Log													
Logging Enabled													
Logging Disabled													
Erase Log													
12:00:00	001/500 												
<p>4. Press . The screen will display Storing... for about 1 second and return to the Editing Menu.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Set Clock</td> <td></td> </tr> <tr> <td>Logging</td> <td></td> </tr> <tr> <td>Factory Setup</td> <td></td> </tr> <tr> <td>Set Power Save</td> <td></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Editing Menu		Set Clock		Logging		Factory Setup		Set Power Save		12:00:00	001/500 
Editing Menu													
Set Clock													
Logging													
Factory Setup													
Set Power Save													
12:00:00	001/500 												

■ FACTORY SETUP

The Factory Setup menu is used in manufacturing of the colorimeter. This menu is not for use by the operator in the field.

■ SETTING POWER SAVE





























The power saving Auto Shutoff feature will turn the meter off when a button has not been pushed for a set amount of time. The default setting is disabled. To change the setting:

<p>1. From the Editing Menu, press  or  to scroll to Set Power Save.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Set Clock</td> <td rowspan="4" style="text-align: center; vertical-align: middle;">   </td> </tr> <tr> <td>Logging</td> </tr> <tr> <td>Factory Setup</td> </tr> <tr> <td>Set Power Save</td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Editing Menu		Set Clock	 	Logging	Factory Setup	Set Power Save	12:00:00 001/500 				
Editing Menu													
Set Clock	 												
Logging													
Factory Setup													
Set Power Save													
12:00:00 001/500 													
<p>2. Press  to select Set Power Save.</p>	<table border="1"> <thead> <tr> <th colspan="2">Set Power Save</th> </tr> </thead> <tbody> <tr> <td>Disable</td> <td style="text-align: center;"> </td> </tr> <tr> <td>5 Minutes</td> <td style="text-align: center;"> </td> </tr> <tr> <td>15 Minutes</td> <td style="text-align: center;"> </td> </tr> <tr> <td>30 Minutes</td> <td style="text-align: center;"> </td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Set Power Save		Disable		5 Minutes		15 Minutes		30 Minutes		12:00:00 001/500 	
Set Power Save													
Disable													
5 Minutes													
15 Minutes													
30 Minutes													
12:00:00 001/500 													
<p>3. Press  or  to scroll to desired function.</p>	<table border="1"> <thead> <tr> <th colspan="2">Set Power Save</th> </tr> </thead> <tbody> <tr> <td>Disable</td> <td style="text-align: center;"> </td> </tr> <tr> <td>5 Minutes</td> <td style="text-align: center;"> </td> </tr> <tr> <td>15 Minutes</td> <td style="text-align: center;"> </td> </tr> <tr> <td>30 Minutes</td> <td style="text-align: center;"> </td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Set Power Save		Disable		5 Minutes		15 Minutes		30 Minutes		12:00:00 001/500 	
Set Power Save													
Disable													
5 Minutes													
15 Minutes													
30 Minutes													
12:00:00 001/500 													
<p>4. Press . The screen will display Storing.... for about 1 second and the meter will return to the Editing Menu.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Set Clock</td> <td rowspan="4" style="text-align: center; vertical-align: middle;">   </td> </tr> <tr> <td>Logging</td> </tr> <tr> <td>Factory Setup</td> </tr> <tr> <td>Set Power Save</td> </tr> <tr> <td colspan="2">12:00:00 001/500 </td> </tr> </tbody> </table>	Editing Menu		Set Clock	 	Logging	Factory Setup	Set Power Save	12:00:00 001/500 				
Editing Menu													
Set Clock	 												
Logging													
Factory Setup													
Set Power Save													
12:00:00 001/500 													

■ SETTING THE BACKLIGHT TIME





The backlight illuminates the display for enhanced viewing. The default setting is 10 seconds. If Button Control is chosen the backlight button on the key pad will act as an on/off switch and the backlight will remain on or off when the meter is being used. When one of the other settings – 10, 20 or 30 seconds – is chosen, the display will be illuminated for the specified amount of time after any button is pressed.




NOTE: The backlight feature uses a significant amount of power. The longer the backlight is on, the more frequently the battery will have to be charged if the USB/Wall Adapter is not being used.

<p>1. From the Editing Menu, press  or  to scroll to Backlight Time.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Logging</td> <td></td> </tr> <tr> <td>Factory Setup</td> <td></td> </tr> <tr> <td>Set Power Save</td> <td></td> </tr> <tr> <td>Set Backlight Time</td> <td></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Editing Menu		Logging		Factory Setup		Set Power Save		Set Backlight Time		12:00:00	001/500 
Editing Menu													
Logging													
Factory Setup													
Set Power Save													
Set Backlight Time													
12:00:00	001/500 												
<p>2. Press ENTER to select Set Backlight Time.</p>	<table border="1"> <thead> <tr> <th colspan="2">Set Backlight Time</th> </tr> </thead> <tbody> <tr> <td>Button Control</td> <td><input type="checkbox"/></td> </tr> <tr> <td>10 seconds</td> <td><input type="checkbox"/></td> </tr> <tr> <td>20 seconds</td> <td><input type="checkbox"/></td> </tr> <tr> <td>30 seconds</td> <td><input type="checkbox"/></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Set Backlight Time		Button Control	<input type="checkbox"/>	10 seconds	<input type="checkbox"/>	20 seconds	<input type="checkbox"/>	30 seconds	<input type="checkbox"/>	12:00:00	001/500 
Set Backlight Time													
Button Control	<input type="checkbox"/>												
10 seconds	<input type="checkbox"/>												
20 seconds	<input type="checkbox"/>												
30 seconds	<input type="checkbox"/>												
12:00:00	001/500 												
<p>3. Press  or  to scroll to desired option.</p>	<table border="1"> <thead> <tr> <th colspan="2">Set Backlight Time</th> </tr> </thead> <tbody> <tr> <td>Button Control</td> <td><input type="checkbox"/></td> </tr> <tr> <td>10 seconds</td> <td><input type="checkbox"/></td> </tr> <tr> <td>20 seconds</td> <td><input checked="" type="checkbox"/></td> </tr> <tr> <td>30 seconds</td> <td><input type="checkbox"/></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Set Backlight Time		Button Control	<input type="checkbox"/>	10 seconds	<input type="checkbox"/>	20 seconds	<input checked="" type="checkbox"/>	30 seconds	<input type="checkbox"/>	12:00:00	001/500 
Set Backlight Time													
Button Control	<input type="checkbox"/>												
10 seconds	<input type="checkbox"/>												
20 seconds	<input checked="" type="checkbox"/>												
30 seconds	<input type="checkbox"/>												
12:00:00	001/500 												
<p>4. Press ENTER. The screen will display Storing... for about 1 second and the meter will return to the Editing Menu.</p>	<table border="1"> <thead> <tr> <th colspan="2">Editing Menu</th> </tr> </thead> <tbody> <tr> <td>Logging</td> <td></td> </tr> <tr> <td>Factory Setup</td> <td></td> </tr> <tr> <td>Set Power Save</td> <td></td> </tr> <tr> <td>Set Backlight Time</td> <td></td> </tr> <tr> <td>12:00:00</td> <td>001/500 </td> </tr> </tbody> </table>	Editing Menu		Logging		Factory Setup		Set Power Save		Set Backlight Time		12:00:00	001/500 
Editing Menu													
Logging													
Factory Setup													
Set Power Save													
Set Backlight Time													
12:00:00	001/500 												




■ SELECTING A LANGUAGE

There are seven languages available: English, Spanish, French, Portuguese, Italian, Chinese, and Japanese.

1. From the Editing Menu , press  or  to scroll to Select Language .	Editing Menu
	<div style="background-color: #333; color: #fff; padding: 5px;"> Factory Setup Set Power Save Set Backlight Time Select Language </div> <div style="text-align: right; margin-top: -10px;"></div>
	12:00:00 001/500 






2. Press  to select Select Language .	Select Language
	<div style="background-color: #333; color: #fff; padding: 5px;"> English Spanish French Portugese </div> <div style="text-align: right; margin-top: -10px;"></div>
	12:00:00 001/500 

3. Press  or  to scroll to desired language.	Select Language
	<div style="background-color: #333; color: #fff; padding: 5px;"> English Spanish French Portugese </div> <div style="text-align: right; margin-top: -10px;"></div>
	12:00:00 001/500 

4. Press  . The screen will display Storing... for about 1 second and the meter will return to the Editing Menu .	Editing Menu
	<div style="background-color: #333; color: #fff; padding: 5px;"> Set Power Save Set Backlight Time Bluetooth Menu Select Language </div> <div style="text-align: right; margin-top: -10px;"></div>
	12:00:00 001/500 


NOTE: If meter unintentionally switches to another language, use the procedure above to reset the meter to the desired language. For example, to reset the meter to English:

Turn meter on.

1. Press  one time. Press .
2. Press  five times. Press .
3. Press .

■ LOOPING MENUS

Long menus, such as All Tests, incorporate a looping feature which allows the user to quickly reach the last choice in the menu from the first choice. In a looping menu the last choices in the menu are above the first choice and scrolling upward moves through the menu in reverse order. Scrolling downward moves through the menu from first choice to last but the menu starts over following the last choice. So all menu choices can be reached by scrolling in either direction. The diagrams below demonstrate a looping menu.

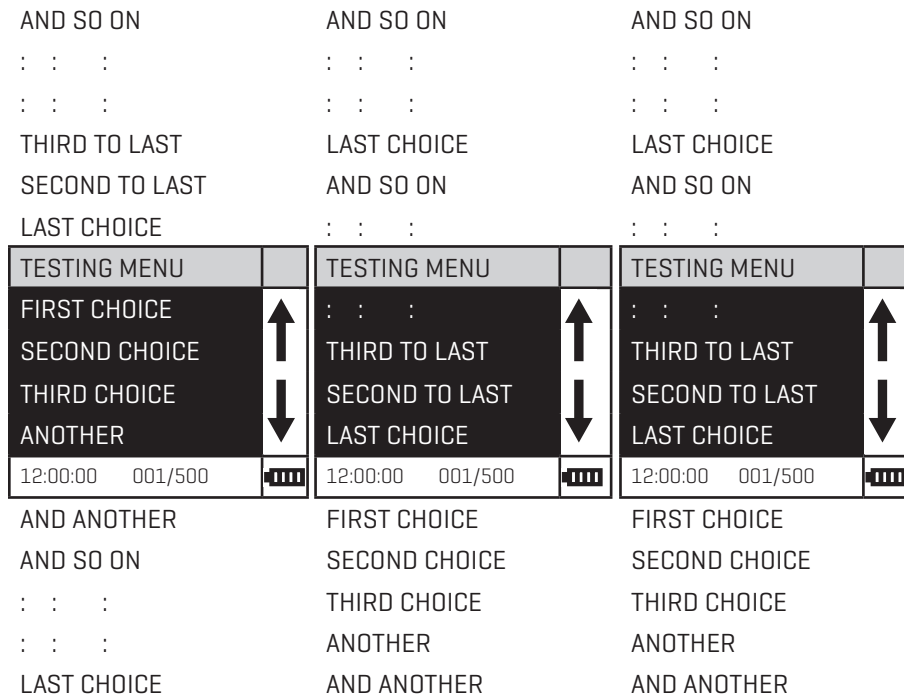

The feature called **Looping Menu** can be turned on and off in the **Editing Menu**. The default setting is ON.

COMPUTER CONNECTION

■ OUTPUT

USB

■ COMPUTER CONNECTION

USB Type A, USB mini B, Order Cable Code 1720.

■ WATERLINK CONNECT

The meter may be interfaced with any Windows-based 64 bit computer by using the LaMotte WaterLink Connect program and a USB cable. The program will store test information and results in a local database, and allow for exporting this data to a comma separated value [CSV] file. The meter will send the following data: time/date stamp, name of test, sample value, sample units, meter name, and location. To download WaterLink go t

BATTERY

■ BATTERY/AC OPERATION

The colorimeter may be operated on battery power using a USB wall adapter or USB computer connection. If using the meter as a bench top unit, use the wall adapter if possible to extend the battery life. The meter will remain on when the USB adapter is used.

To charge the battery with the wall adapter, plug the smaller end of the USB cable [USB mini B connector] into the meter and the larger end of the USB cable [USB type A connector] into the wall adapter. Plug the wall adapter into an AC outlet. Reinsert the rubber USB port plug after charging. To charge the battery from a computer, plug the smaller end of the USB cable [USB mini B connector] into the meter and the larger end of the USB cable [USB Type A connector] into a USB port on a computer. The battery icon will show no bars and flash when the unit first turns on. Then the indicator will indicate the battery status by showing 0, 1, 2, 3, or 4 bars.

It will take 6 hours to fully charge a low battery. The battery icon will flash when the battery is charging. The battery icon will show four bars and stop flashing when it is fully charged. The charging circuit will automatically switch to a float charge when the battery is fully charged. The charger may remain connected. Some computers will NOT supply power to their USB ports during standby operation. The wall adapter will charge the unit continuously.

Storing the meter above ambient room temperature will decrease the battery charge more quickly than storage at room temperature. If the meter does not turn on, it means that the battery is at a very low charge. Charging the battery with the wall adapter in this state may take up to 10 hours. At low temperatures, approaching 0 °C, the battery will charge more slowly. It will not charge at all below 0 °C. The battery icon will show no bars and continuously flash if the battery is getting low but the unit will still operate normally. A "Low Battery" message on the status bar of the display will replace the time

when the battery voltage is too low for proper operation and accuracy may be degraded. A “Shutdown Low Batt” message on the display will appear for a few seconds before the power is switched off when the battery is too low to operate the unit. When the battery icon simultaneously flashes bars 1 and 2 followed by bars 3 and 4, it is an indication that the battery is damaged and technical support should be contacted.

To extend the battery life:

- Shut down the unit with the power switch when not taking measurements or use the power save option to have the unit automatically turn off after 5 minutes.
- Store the unit at a moderate temperature.
- Fully charge the battery before storing the unit for extended periods of time.
- Fully charge the battery at least once per year. Failure to do so may result in a permanently drained battery.
- Limit backlight use. The unit consumes 3X normal power with the backlight on. Set the backlight time option to 10 seconds, or select “Button Control” and keep the backlight off.

BATTERY REPLACEMENT

The lithium ion battery used in this unit should last for many years with normal use. When it no longer powers the unit long enough to meet testing requirements it will need to be replaced. Lithium ion batteries that are properly charged and stored do not usually lose all capacity; they just have less capacity after hundreds of charge cycles. This unit uses a custom battery assembly that is only available from LaMotte Company. Battery replacement must be performed at a LaMotte authorized repair facility. The water resistant housing of this meter should not be opened by the user

MAINTENANCE

■ CLEANING

Clean the exterior housing with a damp, lint-free cloth. Do not allow water to enter the light chamber or any other parts of the meter. To clean the light chamber and optics area, point a can of compressed air into the light chamber and blow the pressurized air into the light chamber. Use a cotton swab dampened with Windex® window cleaner to gently swab the interior of the chamber. Do not use alcohol; it will leave a thin residue over the optics when dry.

■ METER DISPOSAL

Waste Electrical and Electronic Equipment (WEEE)

Natural resources were used in the production of this equipment. This equipment may contain materials that are hazardous to health and the environment. To avoid harm to the environment and natural resources, the use of appropriate take-back systems is recommended. The crossed out wheeled bin symbol on the meter encourages the use of these systems when disposing of this equipment.



Take-back systems will allow the materials to be reused or recycled in a way that will not harm the environment. For more information on approved collection, reuse, and recycling systems contact local or regional waste administration or recycling services.

GENERAL OPERATING INFORMATION

■ OVERVIEW

The meter is a portable, microprocessor controlled, direct reading colorimeter. It has a graphical liquid crystal display and 6 button keypad. These allow the user to select options from the menu driven software, to directly read test results or to review stored results of previous tests in the data logger. The menus can be displayed in seven different languages.

The LaMotte tests are precalibrated for LaMotte reagent systems. The colorimeter displays the result of these tests directly in units of concentration.

The optics feature a colored LED. The LED has a corresponding silicon photoiode with an integrated interference filter. The interference filter selects a narrow band of light from the corresponding LED for the colorimetric measurements.







A USB wall adapter, USB computer connection or lithium battery powers the colorimeter.


GENERAL OPERATING INFORMATION

The operation of the colorimeter is controlled by the menu driven software and user interface. A menu is a list of choices. This allows a selection of various tasks for the colorimeter to perform, such as scan blank and scan sample. The keypad is used to make menu selections that are viewed on the display.

■ THE KEYPAD

The keypad has 6 buttons which are used to perform specific tasks.

	This button will scroll up through a list of menu selections.
	The button is used to select choices in a menu viewed on the display.
	This button controls the backlight on the display.
	This button will scroll down through a list of menu selections.
	This button exits to the previous menu.
	This button turns the meter on or off.

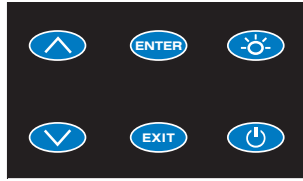


■ THE DISPLAY AND MENU

The display allows menu selections to be viewed and selected. These selections instruct the colorimeter to perform specific tasks. The menus are viewed in the display using two general formats that are followed from one menu to the next. Each menu is a list of choices or selections.

The display has a header line at the top and a footer line at the bottom. The header displays the title of the current menu. The footer line displays the time and the date, the data logger status and the battery status. The menu selection window is in the middle of the display between the header and the footer.


The menu selection window displays information in two general formats. In the first format only menu selections are displayed. Up to 4 lines of menu selections may be displayed. If more selections are available they can be viewed by pressing the arrow buttons to scroll the other menu selections into the menu selection window. Think of the menu selections as a vertical list in the display that moves up or down each time an arrow button is pressed. Some menus in the colorimeter are looping menus. The top and bottom menu choices are connected in a loop. Scrolling down past the bottom of the menu will lead to the top of the menu. Scrolling up past the top of the menu will lead to the bottom of the menu.

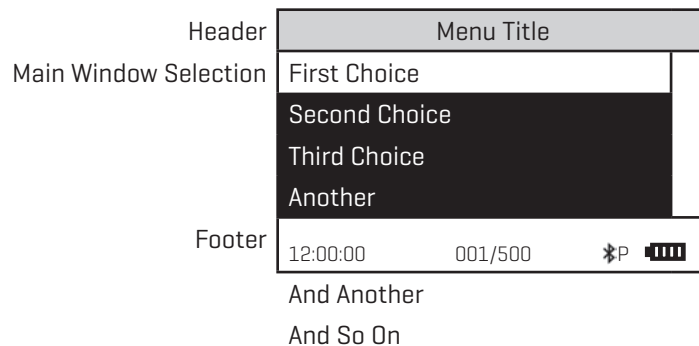


A light bar will indicate the menu choice. As the menu is scrolled through, the light bar will highlight different menu choices. Pressing the button will select the menu choice that is indicated by the light bar.



In the second format the menu choice window takes advantage of the graphical capabilities of the display. Large format graphic information, such as test results or error messages or the LaMotte logo is displayed. The top two lines of the display are used to display information in a large, easy to read

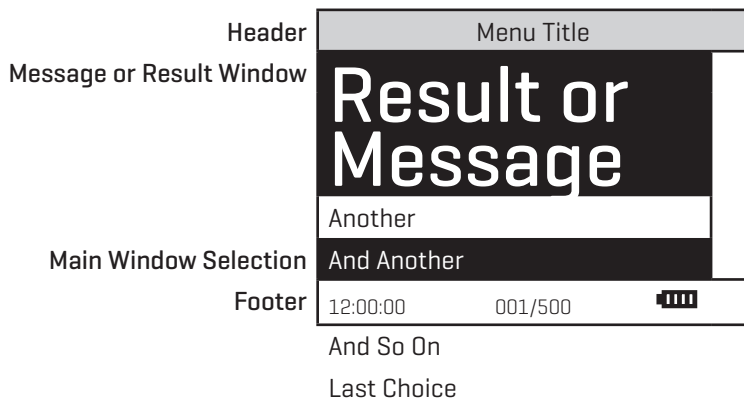format. The menus work in the same way as previously described but two lines of the menu are visible at the bottom of the display.



As described previously, the **EXIT** button allows an exit or escape from the current menu and a return to the previous menu. This allows a rapid exit from an inner menu to the main menu by repeatedly pushing the **EXIT** button. Pushing  at any time will turn the colorimeter off.

The display may show the following messages:

	Battery Status
	More choices are available and can be viewed by scrolling up and/or down through the display.
Header	Identifies the current menu and information on units and reagent systems if applicable.
Footer	In the data logging mode the number of the data point is displayed and the total number of data points in the memory will be shown. The footer also shows current time and battery status

■ TUBES AND CHAMBERS

The colorimeter uses one type of tube [Code 0290] for all test factors.

The handling of the tubes is of utmost importance. Tubes must be clean and free from lint, fingerprints, dried spills and significant scratches, especially the central zone between the bottom and the sample line.

Scratches, fingerprints and water droplets on the tube can cause stray light interference leading to inaccurate results. Tubes that have been scratched in the light zone through excessive use should be discarded and replaced with new ones.

Tubes should always be washed on the inside and outside with mild detergent prior to use to remove dirt or fingerprints. The tubes should be allowed to air-dry in an inverted position to prevent dust from entering the tubes. Dry tubes should be stored with the caps on to prevent contamination.

After a tube has been filled and capped, it should be held by the cap and the outside surface should be wiped with a clean, lint-free absorbent cloth until it is dry and smudge-free. Handling the tube only by the cap will avoid problems from fingerprints. Always set the clean tube aside on a clean surface that will not contaminate the tube. It is imperative that the tubes and light chamber be clean and dry. The outside of the tubes should be dried with a clean, lint-free cloth or disposable wipe before they are placed in the meter chamber.

Tubes should be emptied and cleaned as soon as possible after reading a sample to prevent deposition of particulates on the inside of the tubes.

Variability in the geometry of the glassware and technique is the predominate cause of variability in results. Slight variations in wall thickness and the diameter of the tubes may lead to slight variations in the test results. To eliminate this error the tubes should be placed in the chamber with the same orientation each time.

Chambers which have been scratched through excessive use should be discarded and replaced with a new one.

■ SAMPLE DILUTION TECHNIQUES

If a test result is out of the range of the meter, it must be diluted. The test should then be repeated on the diluted sample. The following table gives quick reference guidelines for dilutions of various proportions.

Amount of Sample	Deionized Water to Bring Final Volume to 10 mL	Multiplication Factor
10 mL	0 mL	1
5 mL	5 mL	2
2.5 mL	7.5 mL	4
1 mL	9 mL	10
0.5 mL	9.5 mL	20

All dilutions are based on a final volume of 10 mL, so several dilutions will require small volumes of the water sample. Graduated pipets should be used for all dilutions. If volumetric glassware is not available, dilutions can

be made with the colorimeter tube. Fill the tube to the 10 mL line with the sample and then transfer it to another container. Add 10 mL volumes of deionized water to the container and mix. Transfer 10 mL of the diluted sample to the colorimeter tube and follow the test procedure. Repeat the dilution and testing procedures until the result falls within the range of the calibration. Multiply the test result by the dilution factor. For example, if 10 mL of the sample water is diluted with three 10 mL volumes of deionized water, the dilution factor is four. The test result of the diluted sample should be multiplied by four.

GENERAL INFORMATION

■ PACKAGING AND DELIVERY

Experienced packaging personnel at LaMotte Company assure adequate protection against normal hazards encountered in transportation of shipments.

After the product leaves LaMotte Company, all responsibility for safe delivery is assured by the transportation company. Damage claims must be filed immediately with the transportation company to receive compensation for damaged goods.

■ GENERAL PRECAUTIONS

READ THE INSTRUCTION MANUAL BEFORE ATTEMPTING TO SET UP OR

OPERATE THE METER. Failure to do so could result in personal injury or damage to the meter. The meter should not be used or stored in a wet or corrosive environment. Care should be taken to prevent water from wet tubes from entering the meter chamber.

NEVER PUT WET TUBES IN THE METER.

■ SAFETY PRECAUTIONS

Read the labels on all LaMotte reagent containers prior to use. Some containers include precautionary notices and first aid information. Certain reagents are considered hazardous substances and are designated with a * in the instruction manual. Safety Data Sheets [SDS] can be found at

Keep equipment and reagent chemicals out of the reach of young children.

Ensure that the protection provided by this equipment is not impaired. Do not install or use this equipment in a manner that is not indicated in this manual.

■ LIMITS OF LIABILITY

Under no circumstances shall LaMotte Company be liable for loss of life, property, profits, or other damages incurred through the use or misuse of its products.

■ SPECIFICATIONS & RANGES

INSTRUMENT TYPE: Colorimeter

Readout	160 x 100 backlit LCD, 20 x 6 line graphical display
Wavelengths	525 nm
Wavelength Accuracy	±2% FS
Readable Resolution	Determined by reagent system
Wavelength Bandwidth	10 nm typical
Photometric Range	-2 to +2 AU
Photometric Precision	± 0.001 AU at 1.0 AU
Photometric Accuracy	±0.005 AU at 1.0 AU
Sample Chamber	Accepts 25 mm diameter flat-bottomed test tubes
Light Sources	1 LEDs
Detectors	1 silicon photodiode
Modes	Pre-programmed tests, absorbance, %T
Pre-Programmed Tests	YES, with automatic wavelength selection
Languages	English, Spanish, French, Portuguese, Italian, Chinese, Japanese
Temperature	Operation: 0-50 °C; Storage: -40-60 °C
Operation Humidity Range	0-90 % RH, non-condensing
USB Port	Mini B
Power Requirements	USB wall adapter, USB computer connection or lithium ion rechargeable battery
Battery	Charge Life: Approximately 380 tests with backlight on to 1000 tests with backlight off. Battery Life: Approximately 500 charges.
Electrical Rating	Rated voltage [5V], Rated power of input current [1.0A] at mini-USB input port
Data Logger	500 test results
Waterproof	IP67 with USB port plug in place
Dimensions [LxWxH]	3.5 x 7.5 x 2.5 inches, 8.84 x 19.05 x 6.35 cm
Weight	13 oz, 362 g [meter only]

■ CE COMPLIANCE

The colorimeter has been independently tested and has earned the European CE Mark of compliance for electromagnetic compatibility and safety. To view certificates of compliance, go to [www.itm.com](#).

■ IP67 Certification

The colorimeter meets IP67 standards for protection against dust and immersion only when the USB port plug is in place. Documentation is available at [www.itm.com](#).

TROUBLESHOOTING


■ ERROR MESSAGES

Over Range

If the message **OVERRANGE** is displayed when scanning a sample, the sample may be over range or under range. If the sample is over range the sample should be diluted and tested again [see Sample Dilution Techniques and Volumetric Measurements, page 27].

If overrange is displayed, press **ENTER** to continue testing on diluted samples.

Note: After pressing **ENTER**, the overrange concentration will be displayed. This concentration is an **approximation only**.

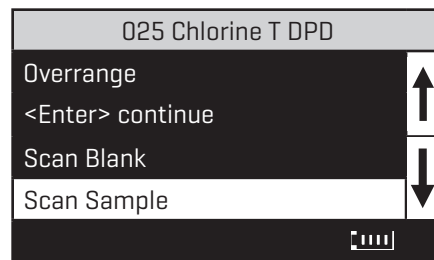


■ CALIBRATION

As with all pre-calibrated meters, it is highly recommended, even if not required by regulations, that the user periodically verify the performance of the meter by running standards with a predetermined concentration. Results outside of specification are an indication that the meter needs to be adjusted. This can be done following the user calibration described on page 24. If the user calibration fails to properly adjust the meter then the meter should be returned to LaMotte Company for recalibration. [See page 30].

■ STRAY LIGHT

The colorimeter should have no problems with stray light. Make sure that the sample compartment lid is always fully closed.

■ TROUBLESHOOTING GUIDE

<i>PROBLEM</i>	<i>REASON</i>	<i>SOLUTION</i>
 Flashing	Low battery. Readings are reliable.	Charge battery or use USB wall/computer adapter.
"Low Battery"	Battery voltage is very low. Readings are not reliable.	Charge battery or use USB wall/computer adapter.
"Shut Down Low Batt" Shut Down	Battery is too low to operate the unit.	Charge battery or use USB wall/computer adapter.
"Overrange"	Sample is outside of acceptable range.	Dilute sample and test again.
Unusually large negative or positive readings when performing calibration	Incorrect standards used to calibrate meter.	Use fresh 0.0 standard in clean tube. Reset meter to factory default settings. Recalibrate meter.



3240-LI-MN-CL 08.19